



切割机

QCUT 350 A

Qcut 350 A 是一台自动湿式砂轮切割机，可切割最大直径为**135毫米**的样品。由于其**低接触切割(lcc)**功能，它可以轻柔地切割各种几何形状、尺寸和材料的不同类型的样品。

该切割机通常用于金相学、材相学、矿物学，以及复合材料（CFC/PMC/MMC）或电子装置（如PCB）的切割。

该机器的大型切割室和不锈钢工作台使操作舒适而灵活。丰富的基于软件的功能，如自动接触检测、多位置进程和垂直切割时的切割轮保护罩平行设置，进一步提高了切割过程的可重复性。



[点击观看视频](#)

优点

- | 横向自动移动
- | 纵向自动移动
- | 电子控制切割力
- | 电子控制移动距离
- | 电子控制的切割力限制装置
- | 储存切割程序
- | 多种样品夹具和夹持系统可供选择
- | 右侧开口可用于放置长的工件

性能指标

电子性能控制

当达到预设的最大功率时，进给速度会降低。这样可以保护机器及其切割部件，防止过载。

精确横向进给

可选配的自动横向进给和序列切割功能适用于精确的平行切割。通过电子控制进行操作。使用研磨杯轮时，电子横向进给也可用作进给轴。

脉冲切割

在切割重型截面和难以切割的材料时，可变脉冲切割系统具有决定性的优势。

多点进程 MPP

利用 "多点进程" (MPP) 功能，您可以在工作台上放置多个试样，并按一个顺序进行切割。通过示教功能，可以简单方便地设置各种切割位置。

自动接触检测

切割过程自动启动后，样品会快速接近。通过接触切割轮，样品将按照预设的切割参数被高效、安全地切割。

新尺寸

加大的切割室和工作台以及平行设置的砂轮护罩提高了所有轴的切割能力和行程。侧门确保了切割室的最佳进出通道，而用于提供长型样品部件的额外开口则提供了更多可能性。

电子控制

自动横切的切割参数可保存在电子控制装置中，之后再加载。手动横切的位置也会显示在显示屏上。所有参数，如进给速度、行程和切割力，均可在自动切割过程中进行更改。

切割机 QCUT 350 A

超凡细节

1. 大尺寸切割试台

- | 轴可以无反弹定位
- | 简洁的凹槽设计，便于 QATM 夹具的安装
- | 用于单独夹具的12毫米T型槽
- | 最大宽 x 高: 500 mm x 300 mm (无Z轴), 除此外最大 400 mm x 300

2. 人体工学控制面板

- | 10,1" 彩色触摸屏
- | 用于手动操作的操纵杆

3. 激光辅助定位

- | 用于切割前工件的精确定位

4. 切割轮锁定

- | 通过锁定机制和简易螺

5. 侧开口

- | 右侧的开口可以用于放置长的工件

6. 坚固的外壳

- | 坚固的粉末涂层铝质结构

7. 平行导向的切割轮罩盖

- | 交叉切割能力的提升




母系统实现简化的切割
轮更换操作

完美的切割 - 由QPREP耗材保证

当使用正确的元素: 知识、保养和正确选择耗材和切割机进行加工时, 切割更可靠、更安全、更高效。


快速、强大、灵活-QTOOLS 夹具

我们创新的夹具用于安全、精确地夹紧工件, 以确保精确的切割和磨削过程。针对特定的应用和工件尺寸, 我们已经开发了各种不同的夹具。

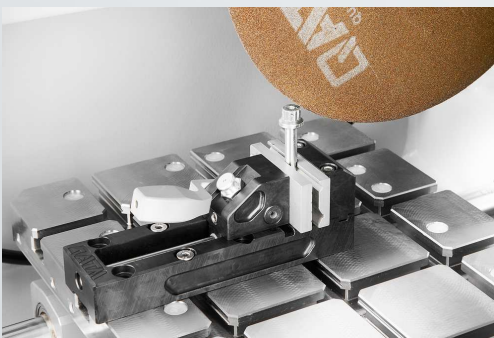

切割机 QCUT 350 A

技术参数

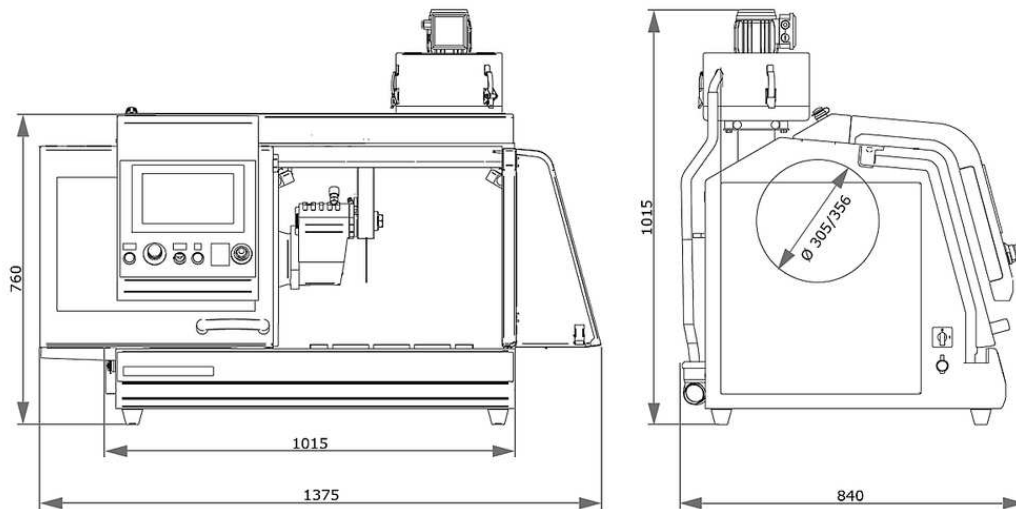

切割轮	up to Ø 356 mm/14"
最大切割能力	Ø 135 mm
切割轮法兰尺寸	Ø 32 mm
自动X轴移动	260 mm
自动Y轴移动	175 mm
Z-轴自动 (可选)	140 mm
Z轴精度	0.1 mm
载物台 (不含Z轴) 宽x深	500 x 300 mm, 不锈钢
载物台 (含Z轴) 宽x深	400 x 300 mm, 不锈钢
T型槽	12 mm
转速	可变转速800-3400转/分
连接电源	9.5 kVA
工作功率 (主载荷)	4 kW S1; 5.5 kW S3
宽 x 高 x 深	1015 x 760 x 810 mm
重量	~200kg(取决于设备)

www.qatm.cn/qcut-350-a

ORDER DATA

BASIC EQUIPMENT

MACHINE (ALL ITEMS REQUIRED)

M1821001  Qcut 350 A (Brillant 250) Base machine

POWER SUPPLY (1 ITEM REQUIRED)

A1821015 Qcut 350 A power supply 400V (3Ph/PE)

A1821016 Qcut 350 A power supply 480V (3Ph/PE)

A1821017 Qcut 350 A power supply 220V (3Ph/PE)

WHEEL GUARD (1 ITEM REQUIRED)

A1821018 Qcut 350 A wheel guard 305mm 12"

A1821019 Qcut 350 A wheel guard 356mm 14"

CROSS FEED (1 ITEM REQUIRED)

A1821020 Qcut 350 A without Z-axis

A1821021 Qcut 350 A manual Z-axis

A1821022 Qcut 350 A automatic Z-axis

RECIRCULATION SYSTEM (1 ITEM REQUIRED)

A5800507  Recirculation cooling System 45L, mobile, with insertion mechanism

A5800508  Recirculation cooling System 45L, mobile, without folding wheels

A6501009  Belt filter 130 - 400V/50Hz (3Ph/PE)

A6501010  Belt filter 130 - 220V/60Hz (3Ph/PE)

Z6501004 Connection set belt filter 130

OPTIONS (TO BE ORDERED WITH MACHINE)

A1821023	Qcut 350 A laser alignment
A1821024	Qcut 350 A centrifugal separator



PRECISION CLAMPING DEVICES

Z1870035	Sample holder L standard Ø 40 mm open
Z1870037	Sample holder L standard Ø 40 mm closed
Z1870039	Sample holder L Flex Ø 40 mm
Z1870043	Sample holder L parallel clamping vice open
Z1870041	Sample holder L parallel clamping vice closed
Z1870045	Drop shape sample holder L Ø 40 mm
Z1870047	Universal clamping vice L
Z1870049	Matrix clamping vice L
Z1870054	Mini vice L
Z1870056	Double parallel vice L
Z1870058	Adaptor with ball joint L
Z1870064	Reduction ring Ø 25 mm
Z1870063	Reduction ring Ø 30 mm
Z1870071	Reduction ring Ø 40 mm
Z1870067	Reduction ring 1"
Z1870066	Reduction ring 1 1/4"
Z1870065	Reduction ring 1 1/2"
Z1870068	Manual chuck Ø 32 mm
Z1870069	Manual chuck Ø 44 mm
Z1870070	Manual chuck Ø 56 mm

VACUUM CLAMPING VICES

Z1870028	Vacuum sample holder
01870288	Vacuum holder (sacrificial plate)
Z1304600	Vacuum pump 230V/50Hz for vacuum sample holder

Z1304634	Vacuum pump 110V/60Hz for vacuum sample holder
----------	--

QTOOL 60

Z2235200	Qtool 60
02235202	Clamping jaw 32 mm - for Qtool 60
Z2235201	Clamping jaws 60 mm - for Qtool 60
02235212	Clamping jaw 32 mm bevelled - for Qtool 60
02235213	Clamping jaw 32 mm cross prism - for Qtool 60
Z2235202	Clamping jaws 60 mm slotted - for Qtool 60
Z2235203	Clamping jaws 60 mm slotted prism - for Qtool 60

QTOOL 80

Z2231203	Qtool 80 - 100mm
02231243	Clamping jaw 40 mm
Z2231215	Clamping jaws 70 mm
02231245	Clamping jaw 40 mm bevelled
02231246	Clamping jaw 40 mm cross prism
Z2231216	Clamping jaws 70 mm slotted
Z2231217	Clamping jaws 70 mm slotted prism

QTOOL 80 VARIO

Z2231200	Qtool 80 Vario - 100 mm
Z2231210	Workpiece stop
02231200	Clamping jaw, 40 mm - for Qtool 80 Vario
Z2231211	Clamping jaws, 70 mm
02231213	Prism jaw, 40 mm
02231229	Bevelled jaw, 40 mm
02231230	Cross prism jaw, 40 mm
Z2231212	Round bar jaws, 70 mm slotted
Z2231213	Prism jaws, 70 mm slotted
Z2231214	Clamping jaws, 70 mm slotted

EASY-CLAMPING PLATES

Z2236030	Easy-clamping plate S
Z2236031	Easy-clamping plate S with rotating unit
Z2236035	Easy-clamping plate M

ACCESSORIES FOR EASY-CLAMPING PLATES

Z2236032	Easy adaption horizontal
Z2236033	Easy adaption vertical
Z2236034	Easy adaption Qtool 40 - 60
Z2236038	Multitower rotating plate
Z2236039	Multitower
Z2236040	Swivel unit Qtool 40 - 60
Z2236036	Easy-multifunctional plate

CLAMPING ACCESSORIES FOR ROTATING, SWIVELING AND SPECIAL APPLICATIONS

Z2236041	Swivel unit Qtool 60 - 80
Z1304643	Screw clamper with Easy-clamping plate M
Z1304636	Expansion clamping device
Z2236009	Vertical clamp single
Z2236010	Vertical clamp double
Z2234010	Clamping system X-Clamp 3858

CROSS VICES

Z1340010	Cross vice 50 - 100
----------	---------------------



ANGLE AND ROTATING VICES

Z1304001	Angle vice 60
----------	---------------

01304008	Prism jaw angle vice 60
----------	-------------------------

ROTATIG DEVICES

Z1870059	Rotation device type RD1
Z1891001	Rotation device type RD2.1

UNIVERSAL VICES AND ACCESSORIES

Z1350025	Clamping rail
01350022	Prism 120°
01350023	Prism slim
01350024	Prism 90°
01350025	Prism 90° high
Z1350008	 Universal vice Piccolo 54
Z1350010	 Suitcase set universal vice Piccolo 54
Z1350012	 Universal vice Piccolo 100
Z1350002	 Universal vice Mono 61
Z1350004	 Suitcase set universal vice Mono 61
Z1350011	 Universal vice Mono 132
Z1350005	Universal vice Duo
Z1350006	Chain vice set
92005796	Longitudinal v-shoe
92005797	Traverse v-shoe
92005799	Longitudinal bridge
92005800	Traverse bridge
92005798	Standard clamping shoe

Z1350007	Height adapter 60mm Piccolo
Z1350003	Height adapter 74mm Mono
95003172	Clamping arm extension Piccolo 54 and Mono 61
Z1304591	Fixing screw M6 x 25
95003169	Fixing screw Piccolo M8 x 25
95003170	Fixing screw Piccolo M10 x 25
95003168	Fixing screw Mono M10 x 32
91001157	T-slot nut - stainless M6 x 8
91001161	T-slot nut - stainless M8 x 10
91001159	T-slot nut - stainless M8 x 12
91001158	T-slot nut - stainless M10 x 12